

LabSwift/LabStart

Changing the DC-power check

Dok-Nr:

006977.00

Blatt 1/2

1 Introduction

This document describes the replacement of the DC Power ribbon cable.

2 Instruction

1. Loosen the cover screws and remove the cover.



2. Disconnect the 6-pin connector and cut the cable.



3. Turn the unit, loosen the marked screws and remove the power check.



4. Insert the new power-check and screw it on.



<p>LabSwift/LabStart</p> <p>Changing the DC-power check</p>	<p>Dok-Nr:</p> <p>006977.00</p>
	<p>Blatt 2/2</p>

5. Thread in ribbon cable and screw down base.



6. Attach the 6-pin socket and crimp it with pliers.



7. Attach and insert the bracket and screw the unit back on.



3 Change process

Vers.	Beschreibung	Datum	Visum
00	Creation of the document	24.03.21	MV/PP